

CradleTM PRO-ISO



CradleTM PRO-ISO

Introduction

Specially designed for CradleTM Pro- Iso , the Esco Shaking CO₂ incubator integrates CelCradleTM technologies. Sleek and reliable Esco CO₂ incubator offers the best temperature and CO₂ concentration control providing superior protection for your cell culture. Shaking function with tide-motion bioreactor helps semi-automated seeding process with up to 98% seeding rate. Optional centrifuge enables to finalize entire cell processing steps inside closed isolator system.

CradleTM Pro-Iso includes single-use, ready-to-use bioreactor called CelCradleTM capable of high density cell culture. CelCradleTM bioreactor is designed based on the concept of bellow-induced intermittent flow of media and air through matrices, where cells resides. Because of its high efficiency in nutrient and oxygen transfer with low shear force, high aeration and foam-free culture environment, one Celcradle bottle (100 ml matrix volume) is capable of producing cell mass comparable to up to 20 roller bottle-850cm².


Applications

- Protein Production
- Virus Production
- Vaccine Research
- Monoclonal Antibody Production
- Cell Bank
- Autologous Cell Therapy: Personal bioreactor
- Allogenic Cell Therapy
- Phase III Vaccine Production Small Scale


Features

CradleTM Pro: Shaking CO₂ incubator + CelcradleTM Bioreactor

- Specially designed incubator for usage in isolator, using Esco Technology, enables evenly heated incubation environment.

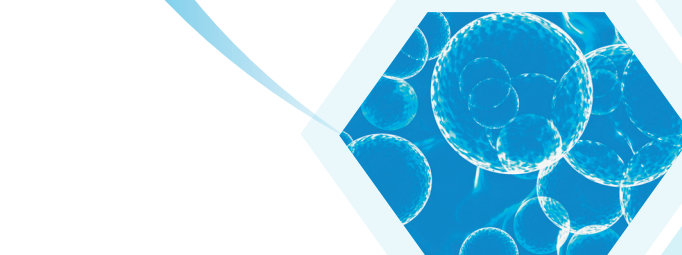


- Able to do pressure test ensures true isolated equipment and environment.
- SteriSafe ULPA filtered system, filtered gas injection lines.
- 0.22 um hydrophobic PTFE-PP filter in the CelcradleTM bottle cap.
- Individual bellow pump and shaking function allow independent use of single CelcradleTM bottle.
- Special treated matrices, BioNOCTM II, enable most anchorage-dependent cells to be cultured.
- Cells entrapped inbetween BioNOCTM II layers simplify media exchange and allow easy harvest and separation of whole cells, cell components and downstream process.
- Optional Stainless Steel 304 exterior is available.

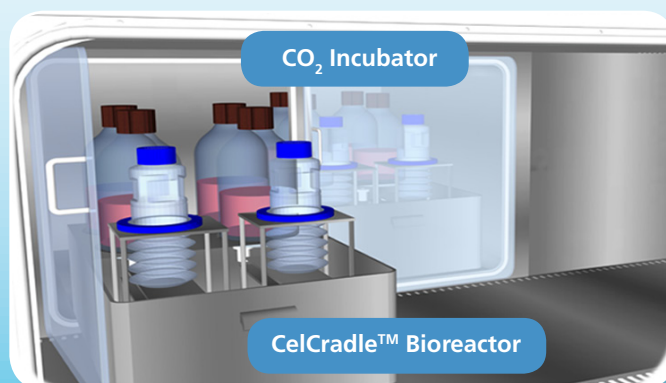

Isolator Specifications



Note: dimensions and components may subject to change depends on further discussion.

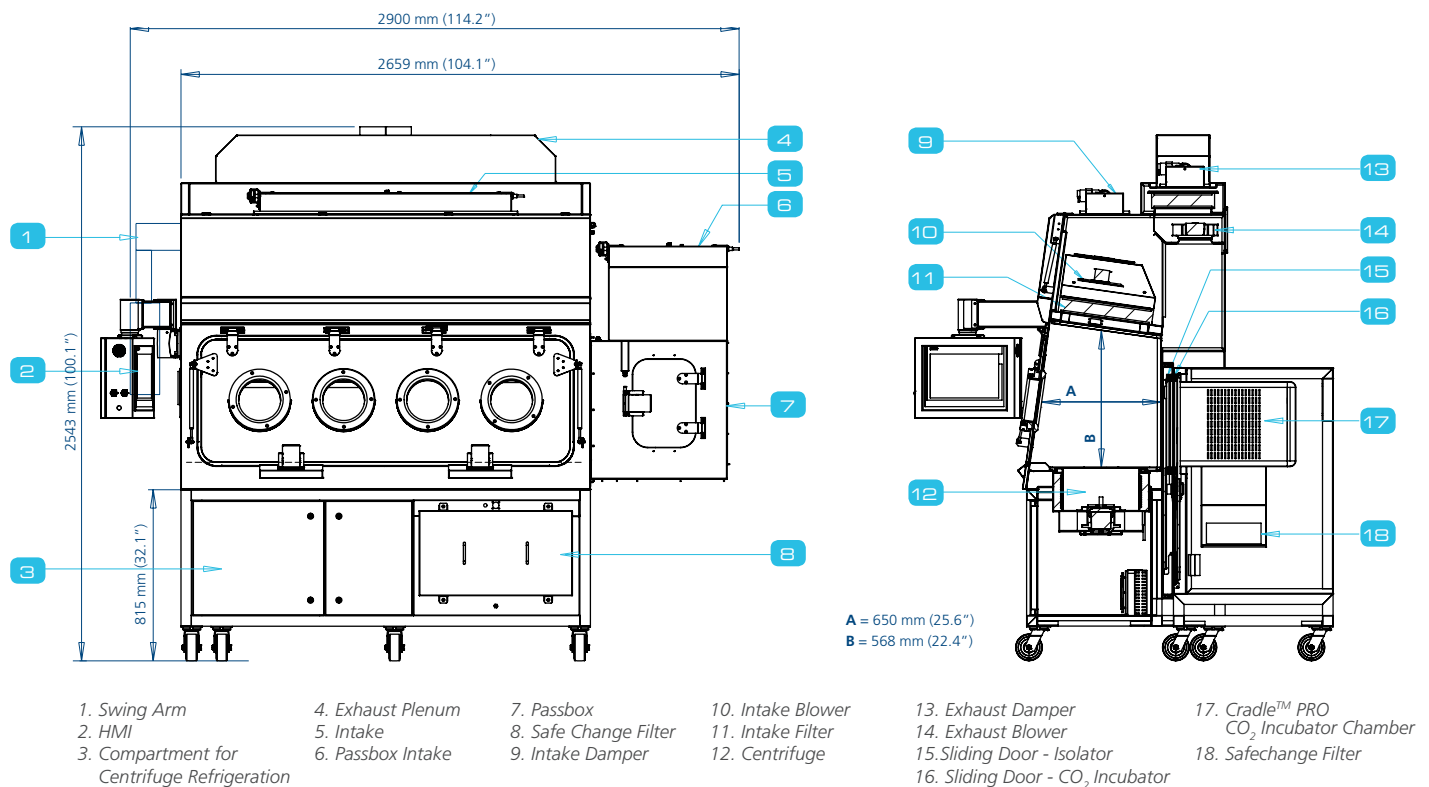

Introduction

- H14 filters with a typical efficiency of 99.999% at 0.1 to 0.3 microns provide superior ISO Class 3 air cleanliness.
- Work zone and pass through chamber are under negative or positive pressure to the room in order to maintain either operator protection or product protection. In case of a breach in the barrier isolation system, there will be either or inward rush in -ve mode or outward air rush in the +ve mode.
- Robust construction and enhanced safety features qualify the Esco Pharmacy Isolator for the most demanding applications.
- Ergonomically angled front and 200mm gloveports improve reach and comfort.
- Esco ISOCIDE™ antimicrobial coating on all painted surfaces minimize contamination. Optional stainless steel 316 L surface possible.
- Cradle™ Pro - Iso is able to be customized depending on client process requirements.

Interior Views

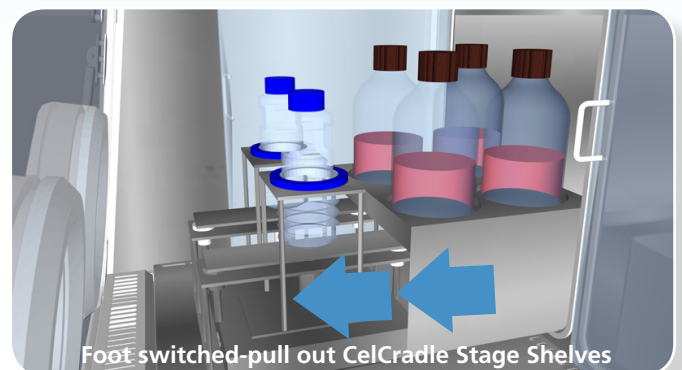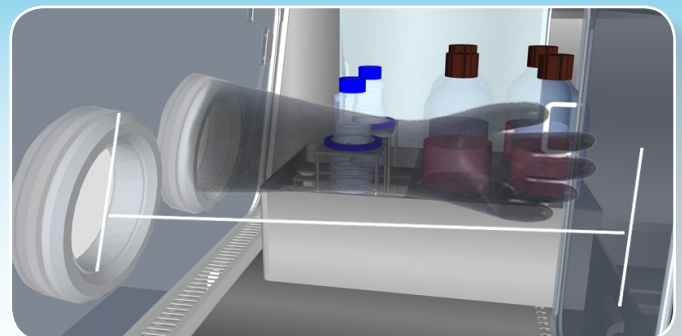

Optional Devices in Esco Isolator

- Microscope to observe cells
- Centrifuge
- Liquid Transfer System
- Benchtop Filling Machine
- Biovap Biodecontamination Hydrogen Peroxide Integration
- Glove Leak Tester